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Thank you,

for purchasing this SCAT Europe product. Please read this operating manual carefully before use and especially observe the safety notices! For service and support, please contact our customer service or your local distributor. For a distributor in your country, please contact us on www.scat-europe.com.

Safety Funnels

Thank you for purchasing our (electrostatic conductive) Safety Funnel. This Funnel has been designed in such a way to comply with the provisions of CENELEC 50404 (Code of practice for the avoidance of hazards due to static electricity) and the German guideline TRGS 727 (formerly TRBS2153) TRGS 727 can be consulted and downloaded on the website www.baua.de. Where to obtain CLC/TR50404 in your country, please request at info@cencenelec.eu

Important notice:

In order to ensure safe operation we strongly recommend to read this operating manual through carefully before putting your funnel into service and to comply with the advices when using it. Please check the chemical resistance of the funnel against the used substances. Also observe the regulations concerning environmental protection and safety in force locally.

Content


Safety Funnel with Hinged Lid and Ball-Valve „ARNOLD“	16
Safety Funnel with Hinged Lid „Marco“	18
Intended use	20
Please notice before any use!	20
Grounding and Antistatic	21
Thermal Properties	23
Personal Protective Equipment.....	23
Maintenance and Care	24
Guarantee and Liability	25

Safety Funnel with Hinged Lid and Ball-Valve „ARNOLD“

Hinged Lid and Tab with Grip

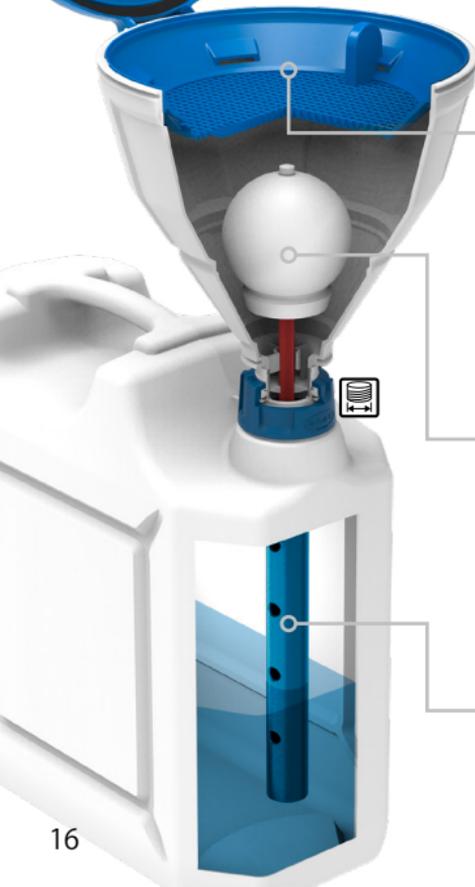All SCAT Europe safety funnels are equipped with the new sealing system that is unique in the world for funnels. Before disposing of liquid waste, open the hinged lid using the tab with grip; the lid remains in its place desired position, so you have both hands free for disposal. After filling the waste canister, close the hinged lid until it clicks into place. The sealing ring closes the funnel cleanly, solvent vapours are retained.

Spray Protection and Dirt Sieve

For collecting mixing magnets or large dirt particles. Laboratory bottles can also be placed here to dry off. The hole size has been optimized, in order to improve the running-off of chemicals. **Both sieve and spray protection can be removed for cleaning.**

Ball-Valve

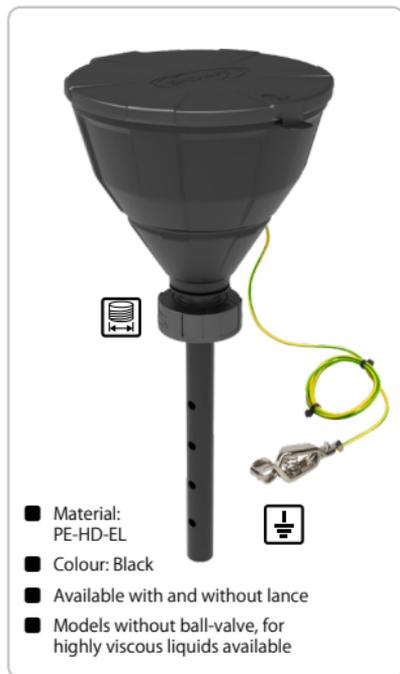
The valve only opens up during filling, and then closes automatically afterwards - for a big „Plus“ regarding safety! To ensure that the ball-valve works properly, please pay attention to the maintenance and care instructions on page 24.

Safety Lance

Ensures for safe and good flowing of sensitive chemicals, without any „stuttering“.

Safety Funnel with Hinged Lid and Ball-Valve „ARNOLD“

EN

The safety funnel „ARNOLD“ is available in various container screw-thread sizes, with and without lance: GL 45, S 50, S 51, B 53, S 55, S 60/61, S 65, S 70/71, B 83, S 90, S 95 and with dual screw thread R2“ BSP/G2“ + 2“ TriSure. In order that we may produce in an environmentally-friendly and economic manner, certain variants are delivered together with a screw-thread adapter. Information for ordering about the funnels, replacement parts as the removable spray protection and dirt sieve, other thread adapters and suitable waste containers you will find at: www.scat-europe.com

Black funnels consist of electrostatic conductive plastic (polyethylene high-density PE-HD) and are produced due to quality management standard ISO 9001. The funnels exhibit the following electrostatic properties: The funnels are electrostatically conductive. The grounding cable and clamp are included in the scope of delivery and already mounted. Please pay attention to the instructions about Grounding and Antistatic on Page 21.

Safety Funnel with Hinged Lid „Marco“

Hinged Lid and Tab with Grip

All SCAT Europe safety funnels are equipped with the new sealing system that is unique in the world for funnels. Before disposing of liquid waste, open the hinged lid using the tab with grip; the lid remains in its place desired position, so you have both hands free for disposal. After filling the waste canister, close the hinged lid until it clicks into place. The sealing ring closes the funnel cleanly, solvent vapours are retained.

Dirt Sieve

For collecting mixing magnets or large dirt particles. Laboratory bottles can also be placed here to dry off. The hole size has been optimized, in order to improve the running-off of chemicals. **The dirt sieve can be removed for cleaning.**

Safety Lance

Ensures for safe and good flowing of sensitive chemicals, without any „stuttering“.

Model with Level Control

Level control integrated in the funnel with float. To ensure that the float works properly, please observe the maintenance and care instructions on page 24.

Safety Funnel with Hinged Lid „Marco“

EN

The safety funnel „Marco“ is available in various container screw-thread sizes, with and without lance: GL 45, S 50, S 51, S 55, S 60/61 and S 65. In order that we may produce in an environmentally-friendly and economic manner, certain variants are delivered together with a screw-thread adapter. Information for ordering about the funnels, replacement parts as the removable dirt sieve, other thread adapters and suitable waste containers you will find at: www.scat-europe.com

Black funnels consist of electrostatic conductive plastic (polyethylene high-density PE-HD) and are produced due to quality management standard ISO 9001. The Funnels exhibit the following electrostatic properties: The funnels are electrostatic conductive. The grounding cable and clamp are included in the scope of delivery and already mounted. Please pay attention to the instructions about Grounding and Antistatic on Page 21.

Intended use

SCAT Europe safety funnels are destined for collecting fluid chemicals, such as solvents, solvent mixtures, water and suspensions. The chemicals may be electrostatic conductive, anti-static or insulating. Any usage beyond this intended use for safety reasons is not allowed.

Funnels with Ball-Valve: By disposing waste with small particles please rinse enough to ensure the proper function of the ball-valve.

Please notice before any use!

When working with the safety funnel, please wear appropriate protective clothing! Always follow the local safety regulations of your company and ask your safety officer in case of any questions!

Please make sure that the safety funnel is installed tightly and free of congestions. When using a safety funnel with ball valve, the ball has to be movable and smooth-running. If the mechanical parts are blocked by dried or viscous fluids, please clean the safety funnel well before using (see „Maintenance and Care“).

Please check the chemical resistance of the funnel material (PE-HD) against the used chemicals. Due to the variety and different constitution of the chemicals available, we do not assume any guarantee or liability regarding chemical resistance. The applicability has to be checked by the operator in each case.

More information on chemical resistance can be provided by the manufacturer of the respective chemicals and can be found in the download area of www.scat-europe.com.

Use the safety funnel only in case of very good chemical resistance. „Very good“ means that the PE-HD material may not show any damage after 30 days of exposure to the chemical.

Do not use the safety funnel in case of insufficient chemical resistance of the PE-HD material. This may cause damage and loss of the electrostatic conductive characteristics.

Grounding and Antistatic

When working with highly flammable fluids with low flashpoint, any device made of electrostatic conductive materials must be grounded safely. Any electrostatic conductive safety funnels made of electrostatic conductive PE-HD (black) must be grounded actively during operation. For grounding, use only the provided grounding cable.

Warning: The installation has to be executed by specialized people only! Please notice the internal safety advises of your company.

The safety funnel is deemed to be grounded if the resistance to ground is max. 10^6 Ohm at a voltage of 500V. If the grounding has been carried out correctly, no electrostatic charges are due when handling electroconductive or antistatic fluids.

When operating with suspensions and insulating fluids, an accumulation of electrostatic charge is possible. This accumulation develops autonomously, independent from the safety funnel. Accumulation on the safety funnel itself may not occur if the grounding has been carried out correctly.

The following factors can affect the electrostatic characteristics of electrostatic conductive safety funnels negatively:

- **The fluid or suspension is sticky and/or insulating.**
This may create a insulating coating at the inner surface of the safety funnel.
- **Diffusion of an insulating fluid** (typically non-conductive solvents) through the plastic material. Over a longer period of time, this may cause loss of the safety funnel's electrostatic conductive characteristics.

Important notice: Electrostatic characteristics of the PE-HD material may change over a longer period of time.

The resistance to the ground has to be checked at least once a year. The resistance to ground has to be max. 10^6 Ohm at a voltage of 500V. In case of deterioration, the safety funnel may no longer be regarded as antistatic and has to be replaced.

Please always follow the local safety regulations of your company and ask your safety officer in case of any questions!

Thermal Properties

The permanent operating temperature of PE without mechanical strain is $-50\text{ }^{\circ}\text{C}$ to $80\text{ }^{\circ}\text{C}$. To avoid the funnel from deformations you should not fill in liquids with a temperature higher than $60\text{ }^{\circ}\text{C}$.

Personal Protective Equipment

Always wear appropriate protective clothing to protect yourself against spillings and vapours during operation.

During maintenance and cleaning of the safety funnel, always wear protective gloves and glasses!

Always follow the local safety regulations of your company and ask your safety officer in case of any questions!

Maintenance and Care

When working with the safety funnel and chemicals, always wear protective clothing and follow the local safety regulations of your company! Ask your safety officer in case of any questions!

After each use, the safety funnel has to be cleaned and purged with an appropriate solvent to avoid congestions or blocking of the ball-valve. In-depth cleaning avoids non-conductive coatings or diffusion that may affect the electrostatic conductive characteristics of the material surface and the anti-static performance of the safety funnel.

Please check the movability of the ball-valve regularly.

Always purge the funnel after using sticky or viscous fluids.

Remove the sieve and the spray protection, and clean them separately.

Purge funnels regularly, e.g. when changing containers, or earlier where necessary. Suitable solvents for purging are water, acetonitrile or isopropyl alcohol. Clean especially movable parts regularly and check them for error-free operation, e.g. the floater when using funnels with level control. This makes sure that there will be no jamming and crystallization caused by drying due to the variety of possible waste liquids.

Please always use wet cloth for external cleaning.

Dry rubbing may cause static charge and sparking when components are not grounded sufficiently.

Guarantee and Liability

Please do not make any technical changes or modifications. If the safety funnel is technically modified by the operator, SCAT Europe GmbH does not assume any guarantee or liability.

If you need a technically modified product for a special application, please contact us. Our product specialists will help you in developing a customized solution.

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The latest version of User Manual relating to your product is always available for you to download. Please visit www.scatt-europe.com or scan this QR-Code using the camera of your Smartphone/Tablet.

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